

AllergyScreen™/AlleisaScreen™ CIRCULAR LETTRE 6


Software-updates for your connect cameras

Software on your PC	You want to update to this software versions	Is it possible to start the installation without a deinstallation of the former version?	structure of database
490	520	yes	old structure
490	530	yes	old structure
490	540	not possible- installation will be interrupted	---
490	540 update	yes	old structure
520	530	yes	new structure
520	540	not possible- installation will be interrupted	---
520	540 update	yes	new structure
530	540	not possible- installation will be interrupted	---
530	540 update	yes	new structure

The new structure of the database was necessary to make it compatible with Windows Vista.

Essential ServicePacks for your PCs to guarantee a correct function of the software

for Windows 2000	Service Pack 4
for Windows XP	Service Pack 3
for Windows Vista	Service Pack 1

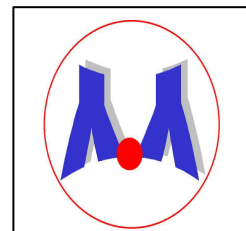

AllergyScreen™/AlleisaScreen™ CIRCULAR LETTRE 6

Reasons for a new software version and the new features

Change in the software versions...	What was the reason for the new version? / What changed?
... from 380 to 437	Preparation for the Logitech 4000 (380 was used for Logitech 3000, 437 was used for Logitech 4000)
... from 437 to 490	Preparation for the Logitech connect (437 was used for Logitech 4000, 490 was the first version, that was used for Logitech connect)
... from 490 to 520	520 is compatible for Windows Vista, Korean language is now available, creating ASCII-files of the day/week is possible with a SINGLE click, gives automatically all rights to new users (doesn't have to be admitted manually), it's possible to print 3 measurements on a sheet, a new designed printout, the clipboard-function for excel-use is available
... from 520 to 530	The software driver was updated
... from 530 to 540	540 was created for a special adaption to the new cameras, greek language is now available, it's possible to delete the headline and footline of the printout to use your own company stationery